



TS-1001	How to Remove the CAB A2+ Printer	Document Owner: Service Department	Effective Date: AUG-23-2022	Page 1 of 5
----------------	--	--	---------------------------------------	--------------------

Purpose

This document provides the steps required to remove and/or replace the Cab A2+ printer on the Sci-Print VX2/VXL.

Intended Use

This document should be used when removing the CAB A2+ printer, and replacing it with the SQUIX2 printer.

Required Material and Equipment

N/A

Intended Audience

Sci-Print VX2/VXL or SOLO Superuser with guidance from a Scinomix Service Technician.

Prerequisites

N/A

Procedure

1. Power down the unit or press to engage the E-Stop.
2. Turn off the air to the unit.
3. Remove the tray by pressing the metal release clip on the tray airline, and gently lift the tray out of the unit.
4. Remove the stainless steel label tensioner by removing the only screw.

TS-1001	How to Remove the CAB A2+ Printer	Document Owner: Service Department	Effective Date: AUG-23-2022	Page 2 of 5
----------------	--	--	---------------------------------------	--------------------

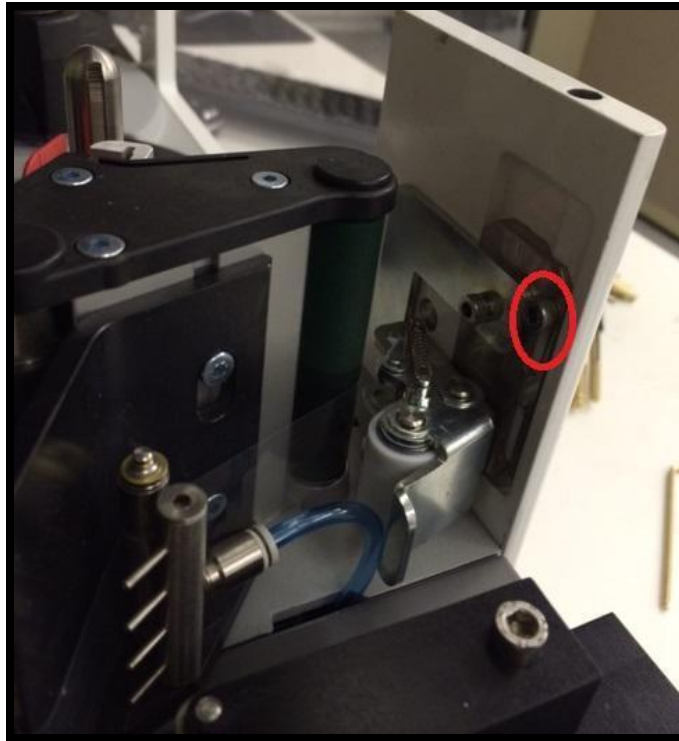


Figure 1.

5. Remove the four screws holding the printer to the base.

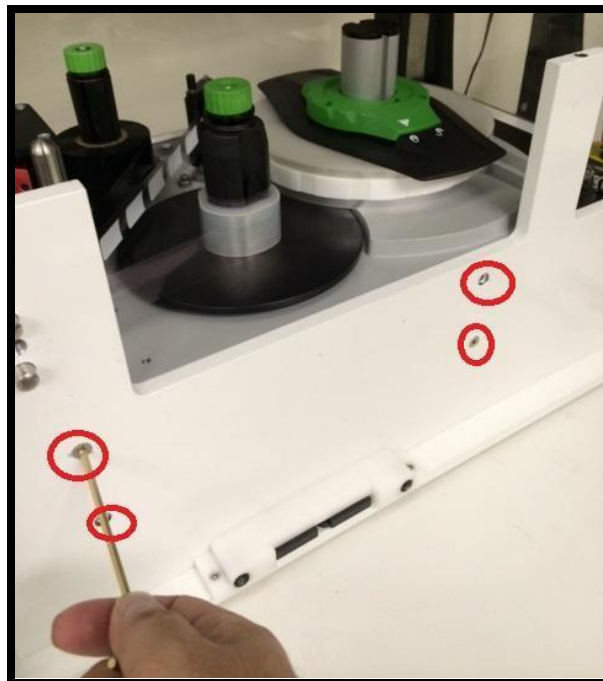


Figure 2.

TS-1001	How to Remove the CAB A2+ Printer	Document Owner: Service Department	Effective Date: AUG-23-2022	Page 3 of 5
----------------	--	--	---------------------------------------	--------------------

6. Turn the run knob on the roller stage to pull the stage out.



Figure 3.

7. Pull the stage out and disconnect the peel blow airline and barcode reader (if equipped).

TS-1001	How to Remove the CAB A2+ Printer	Document Owner: Service Department	Effective Date: AUG-23-2022	Page 4 of 5
----------------	--	--	---------------------------------------	--------------------



Figure 4.

8. Slightly pull the printer out toward the roller stage and unplug the power and Comm line from the back.

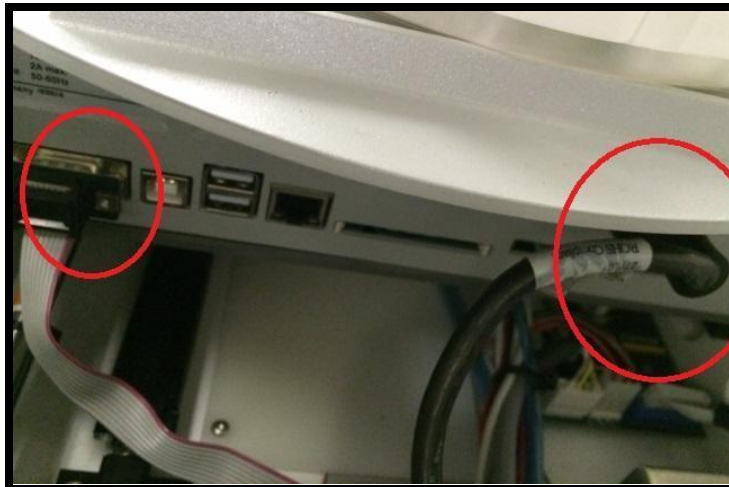


Figure 5.

9. The printer is now free to remove from the unit.
10. Replace with a new printer by following steps in the reverse order.

If you need additional assistance, please call our service department at 314-298-9800 or email service@scinomix.com.



TS-1001	How to Remove the CAB A2+ Printer	Document Owner: Service Department	Effective Date: AUG-23-2022	Page 5 of 5
----------------	--	--	---------------------------------------	--------------------

Definitions

N/A

Revision History

Version:	Change:	Effective Date:	Approved by:
A	NEW	APR-03-2020	B. Petti
B	Process updated to remove A2+ and replace with SQUIX printer	AUG-23-2022	Service