

TS-1005	Issues with VX2-VXL Pneumatic Gripper Crushing Tubes	Document Owner: Service Department	Effective Date: APR-13-2020	Page 1 of 2
----------------	---	--	---------------------------------------	--------------------

Purpose

This guide provides the steps required to correct issues with the VX2/VXL gripper crushing tubes.

Required Material and Equipment

N/A

Intended Audience

Scinomix End user

Prerequisites

Pneumatic grippers are crushing tubes

Procedure

1. Ensure the correct grip finger is being used for the current configuration.
2. Using the stainless steel grip pressure adjustment flow controller, turn the knob counter-clockwise to lessen the pressure.



Figure 1.



TS-1005	Issues with VX2-VXL Pneumatic Gripper Crushing Tubes	Document Owner: Service Department	Effective Date: APR-13-2020	Page 2 of 2
----------------	---	--	---------------------------------------	--------------------

Note: For non-capped tubes, the pressure on the gauge should read no higher than 0.3 MPa. For capped tubes this value should be 0.5 MPa.

If you need additional assistance, please call our service department at 314-298-9800 or email service@scinomix.com.

Definitions

N/A

Revision History

Version:	Change:	Effective Date:	Approved by:
A	NEW	APR-13-2020	B. Petti