

| | | | | |
|----------------|---|--|---------------------------------------|--------------------|
| TS-1014 | How to Replace the CAB A2 Print Head | Document Owner: Service Department | Effective Date: APR-13-2020 | Page 1 of 3 |
|----------------|---|--|---------------------------------------|--------------------|

Purpose

This document provides the steps required to replace the CAB A2 Print Head on the VX2/VXL.

Required Material and Equipment

- 2.5mm allen wrench

Intended Audience

Certified Scinomix Service Technician

Prerequisites

N/A

Procedure

1. Turn off the unit.
2. Open the print head.
3. Remove labels and ribbon from the travel path.
4. Close the print head.
5. Using a 2.5mm allen wrench unscrew the print head from the bracket, then open the print head



Figure 1.

6. Gently pull the print head forward.

| | | | | |
|----------------|---|--|---------------------------------------|--------------------|
| TS-1014 | How to Replace the CAB A2 Print Head | Document Owner: Service Department | Effective Date: APR-13-2020 | Page 2 of 3 |
|----------------|---|--|---------------------------------------|--------------------|

Note: The wires connecting the head can get stuck, so pull on them lightly to free them from behind the bracket.



Figure 2.

7. Unplug the white cable then the black cable, and the print head is free.



Figure 3.

8. Plug in the new print head with the black cable, then the white.
9. Reseat the print head into the print head bracket.
10. Close the print head as much as needed to screw in the bracket screw.



| | | | | |
|----------------|---|--|---------------------------------------|--------------------|
| TS-1014 | How to Replace the CAB A2 Print Head | Document Owner: Service Department | Effective Date: APR-13-2020 | Page 3 of 3 |
|----------------|---|--|---------------------------------------|--------------------|

11. Screw the bracket to the print head using the previously removed screw.
12. Replace labels and ribbon to travel path.
13. Close the print head completely.
14. Turn the unit back on and test.

If you need additional assistance, please call our service department at 314-298-9800 or email service@scinomix.com.

Definitions

N/A

Revision History

| Version: | Change: | Effective Date: | Approved by: |
|-----------------|----------------|------------------------|---------------------|
| A | NEW | APR-13-2020 | B. Petti |