

TS-1050	Removal and Installation of the SQUIX Printer	Document Owner: Service Department	Effective Date: AUG-23-2022	Page 1 of 4
----------------	--	--	---------------------------------------	--------------------

Purpose

This document provides the steps required to remove and install a CAB SQUIX printer.

Intended Use

When a CAB SQUIX needs to be removed and/or replaced.

Required Material and Equipment

- Cab Squix Printer
- 2.5 mm allen wrench

Intended Audience

Superuser

Prerequisites

N/A

Procedure

1. Turn off the instrument.
2. Disconnect the rack lock airline by depressing the silver tab on the quick disconnect air valve while pulling up on the airline connector.



Figure 1.

TS-1050	Removal and Installation of the SQUIX Printer	Document Owner: Service Department	Effective Date: AUG-23-2022	Page 2 of 4
----------------	--	--	---------------------------------------	--------------------

3. Locate and remove label tension assembly spring from the shoulder screw located on the back of the unit base.



Figure 2.

4. Locate and remove the four flathead screws on the front of the base unit with a 2.5 mm Allen wrench.

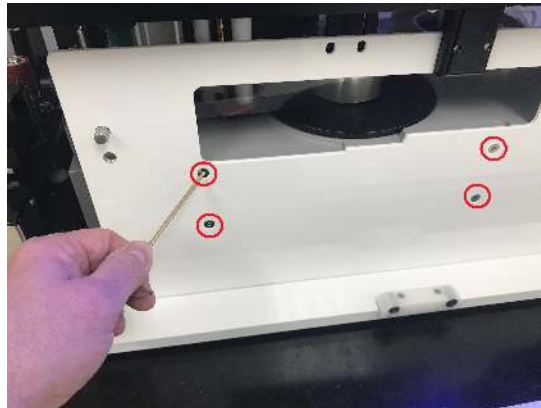


Figure 3.

TS-1050	Removal and Installation of the SQUIX Printer	Document Owner: Service Department	Effective Date: AUG-23-2022	Page 3 of 4
----------------	--	--	---------------------------------------	--------------------

5. Locate and remove the peel blow airline by pushing the small white retaining clip towards the air valve while pulling the airline in the opposite direction.



Figure 4.

6. Turn the run knob ¼ turn to release the roller assembly and pull out the roller to the fully extended position.

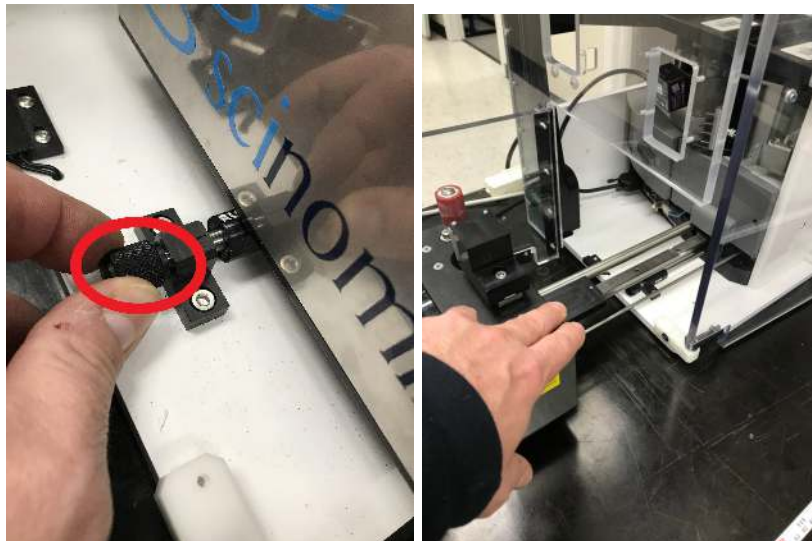


Figure 5.

TS-1050	Removal and Installation of the SQUIX Printer	Document Owner: Service Department	Effective Date: AUG-23-2022	Page 4 of 4
----------------	--	--	---------------------------------------	--------------------

- Slide the printer to the left 4 or 5 inches and lift up on the right side until you can reach in and unplug the power cord and ribbon communication cable.



Figure 6.

- Remove the printer by slowly pulling up making sure not to collide with any other components of the instrument.
- Replace the instrument in the reverse order making sure the Print Power button is in the on position before sliding it into place.

If you need additional assistance, please call our service department at 314-298-9800 or email service@scinomix.com.

Definitions

N/A

Revision History

Version:	Change:	Effective Date:	Approved by:
A	NEW	APR-03-2020	B. Petti
B	Content review <i>Intended Use</i> added	AUG-23-2022	Service