

<b>TS-1053</b>	<b>Grip Finger Jacket Installation for Cryotubes</b>	<b>Document Owner:</b> Service Department	<b>Effective Date:</b> May-05-2020	<b>Page 1 of 2</b>
----------------	--	--	---------------------------------------	--------------------

## Purpose

This document provides the steps required to install new grip finger jackets for cryotubes.

## Required Material and Equipment

N/A

## Intended Audience

Field Service Representative from Scinomix or approved vendors.

## Prerequisites

N/A

## Procedure

1. Push the new grip finger jackets onto the gripper fingers used for cryotubes.
2. Make sure the top of the jacket is against the gripper, to ensure the grip jackets are the same length on the finger.
3. Teach the new position using the **RACK DEFINITION** software application so that the non-jacketed part of the finger is gripping the plastic lip of tube, and the jacketed part of the finger is above the lip of the tube.

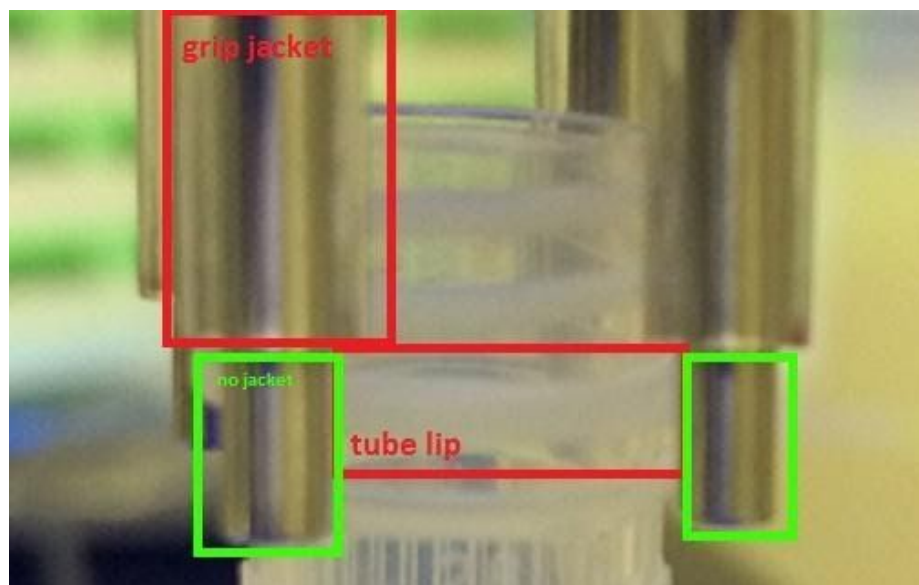


Figure 1.



<b>TS-1053</b>	<b>Grip Finger Jacket Installation for Cryotubes</b>	<b>Document Owner:</b> Service Department	<b>Effective Date:</b> May-05-2020	<b>Page 2 of 2</b>
----------------	--	--	---------------------------------------	--------------------

## Definitions

N/A

## Revision History

<b>Version:</b>	<b>Change:</b>	<b>Effective Date:</b>	<b>Approved by:</b>
A	NEW	MAY-05-2020	B. Petty