

TS-1062	MP2 Label Images	Document Owner: Service Department	Effective Date: MAY-13-2021	Page 1 of 1
----------------	-----------------------------	--	---------------------------------------	--------------------

Purpose

The standard Sci-Print MP2 configuration does not allow printing images on plate labels. This guide offers the steps needed to change this setting by loading a *crc.dll* file which allows images to be printed.

Required Material and Equipment

Sci-Print MP2
crc.dll file from Scinomix

Intended Audience

Scinomix Sci-Print MP2 Users

Prerequisites

N/A

Procedure

1. Download the *crc.dll* or save it to a USB.
2. Right-click on the software icon and select *Open file location*. This shows the location of the install folder.
3. Copy the file from the USB stick (or the download location) and paste it into the install folder (or API folder if integrated).

If you need additional assistance, please call our service department at 314-298-9800 or email service@scinomix.com.

Definitions

N/A

Revision History

Version:	Change:	Effective Date:	Approved by:
A	NEW	MAY-13-2021	J. King