

TS-1068	Updating the Firmware on Squix Printers	Document Owner: Service Team	Effective Date: AUG-24-2022	Page 1 of 4
----------------	--	--	---------------------------------------	--------------------

Purpose

This document provides the steps necessary to update the firmware version on CAB Squix printers.

Intended Use

When new firmware versions are released, requiring an update to the SQUIX printer.

Required Material and Equipment

USB stick with the firmware update

Intended Audience

Certified Scinomix Technician

Prerequisites

For this procedure you need to have the Squix printer powered on, (either on a VX or MP2 instrument, or on a bench). If you use the USB port that is accessible with the printer in the instrument, the USB stick needs to have a slender housing to clear the CAB deck.

Procedure

1. Remove the black USB cover if using the top USB port.



Figure 1.

TS-1068	Updating the Firmware on Squix Printers	Document Owner: Service Team	Effective Date: AUG-24-2022	Page 2 of 4
---------	---	---------------------------------	--------------------------------	-------------

2. Insert the USB stick into either USB port.

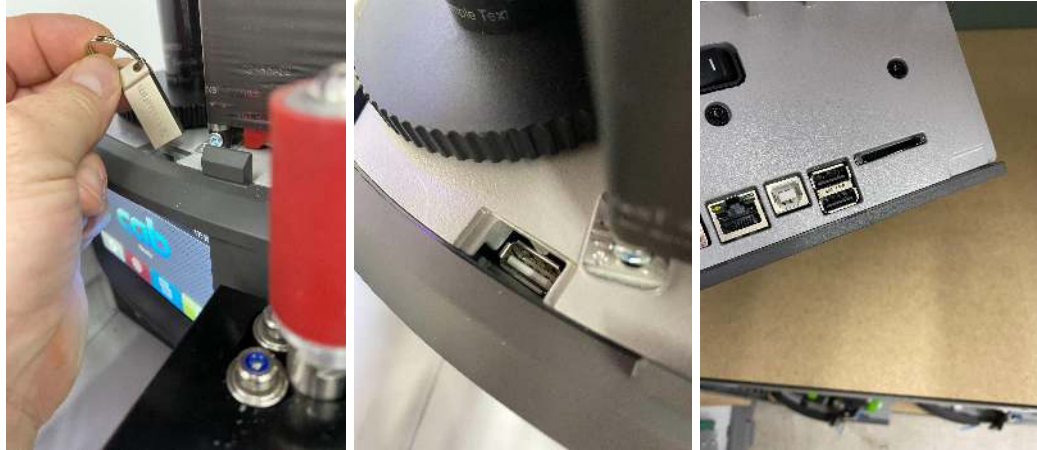


Figure 2.

3. Open the **Settings** tab on the Squix printer LCD screen.
4. Press to open **Extras**.



Figure 3.

TS-1068	Updating the Firmware on Squix Printers	Document Owner: Service Team	Effective Date: AUG-24-2022	Page 3 of 4
---------	---	---------------------------------	--------------------------------	-------------

5. Press **Firmware Update**.



Figure 4.

6. Select and highlight the desired firmware version from the USB stick.



Figure 5.

7. Press the green checkmark in the bottom left of the screen.



Figure 6.

8. The firmware update runs and typically completes within a minute.

9. When complete, open the Home screen and press **Settings** to view the information and verify the firmware version is updated.



TS-1068	Updating the Firmware on Squix Printers	Document Owner: Service Team	Effective Date: AUG-24-2022	Page 4 of 4
----------------	--	--	---------------------------------------	--------------------

If you need additional assistance, please call our Service Department at 314-298-9800 or email service@scinomix.com.

Definitions

N/A

Revision History

Version:	Change:	Effective Date:	Approved by:
A	NEW	JUL-13-2021	M. Tapp
B	Content review Added <i>Intended Use</i>	AUG-24-2022	Service