

<b>TS-1088</b>	<b>Adjusting the Printer Heat Setting (A2+ of VX2/VXL/Solo)</b>	<b>Document Owner:</b> Service Department	<b>Effective Date:</b> AUG-22-2022	<b>Page 1 of 4</b>
----------------	---	--	---------------------------------------	--------------------

## Purpose

This document provides the steps required to adjust the heat setting on the Cab A2+ printer.

## Intended Use

If a different heat setting is needed on the A2+ printer.

## Required Material and Equipment

N/A

## Intended Audience

End users

## Prerequisites

N/A

## Procedure

1. Press the button labeled MENU (this is also the down arrow).



Figure 1.

2. Scroll to SETUP using the button labeled FEED (up arrow).



Figure 2.

<b>TS-1088</b>	<b>Adjusting the Printer Heat Setting (A2+ of VX2/VXL/Solo)</b>	<b>Document Owner:</b> Service Department	<b>Effective Date:</b> AUG-22-2022	<b>Page 2 of 4</b>
----------------	---	--	---------------------------------------	--------------------

3. Press CANCEL (right arrow) to enter the Setup menu.



Figure 3.

4. Scroll to PRINT PARAMS using FEED (up arrow).

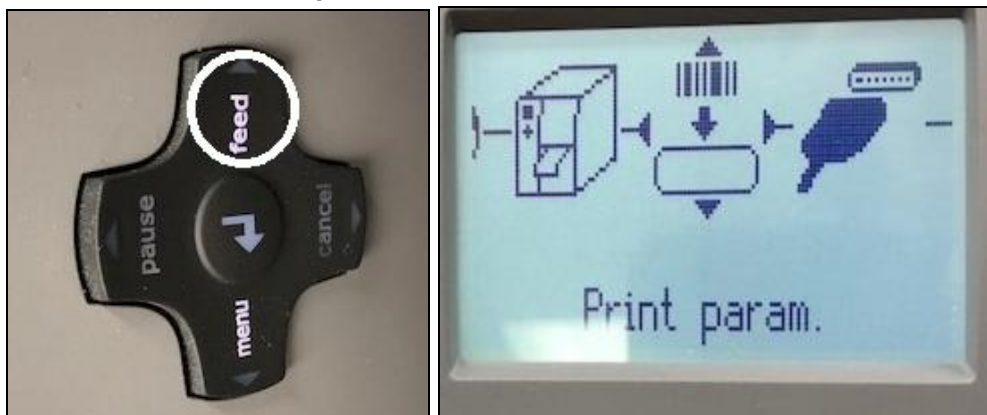


Figure 4.

5. Press CANCEL (right arrow) to enter the Print parameters menu.



Figure 5.

<b>TS-1088</b>	<b>Adjusting the Printer Heat Setting (A2+ of VX2/VXL/Solo)</b>	<b>Document Owner:</b> Service Department	<b>Effective Date:</b> AUG-22-2022	<b>Page 3 of 4</b>
----------------	---	--	---------------------------------------	--------------------

6. The Heat level screen opens.

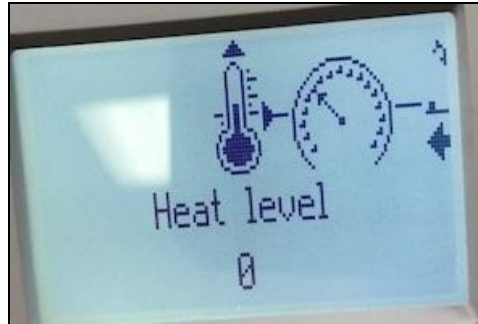


Figure 6.

7. Press the Center button to enter the heat adjustment scale.



Figure 7.

8. Press + until the setting is up to 10.



Figure 8.

9. Press the Center button again to save this setting.



<b>TS-1088</b>	<b>Adjusting the Printer Heat Setting (A2+ of VX2/VXL/Solo)</b>	<b>Document Owner:</b> Service Department	<b>Effective Date:</b> AUG-22-2022	<b>Page 4 of 4</b>
----------------	---	--	---------------------------------------	--------------------

10. HOLD the Center button down to return to a Ready screen.

*If you need additional assistance, please call our service department at 314-298-9800 or email [service@scinomix.com](mailto:service@scinomix.com).*

## Definitions

N/A

## Revision History

<b>Version:</b>	<b>Change:</b>	<b>Effective Date:</b>	<b>Approved by:</b>
A	New	AUG-22-2022	Service